



FAB^{ion}[®]

CELL SELECTION REINVENTED

> 90% PURITY
FROM
WHOLE BLOOD




CELL SELECTION FROM WHOLE

FAB^{ion}®

FAB-BASED

IMMUNE AFFINITY CHROMATOGRAPHY

- › fast
- › from whole blood
- › high yields
- › high purity
- › viability: >99%
- › magnetic bead free
- › fully automatic
- › embedded system
- › label-free authentic target cells
- › no centrifugation
- › web-based programming
- › user independent

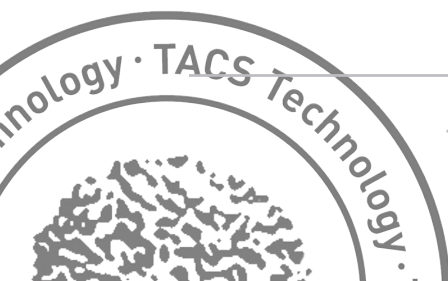

THE SELECTION PROCESS OUTCOME: LABEL FREE, PURI





WHOLE BLOOD MAGNETIC BEAD FREE

CELL SELECTION FROM WHOLE BLOOD

The FABian® cell selection process employs IBA's unique traceless affinity cell selection (TACS) technology for the automatic selection of specialized cell populations including T & B cells and other cell types.

The novel cell selection technology utilizes non-magnetic and fully reversible Strep-tagged Fab fragments targeting specific cell surface markers (e.g. CD81, CD3, CD4, CD8, CD14 and CD19).

FABian® quantitatively selects the cells of interest in high yields with high efficiency and purity from a suspension such as whole blood, buffy coat and other single-cell tissue suspensions. Due to the mild procedure to gain label-free target cells, serial positive cell selections are possible.

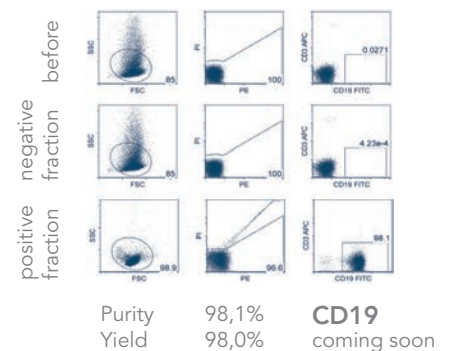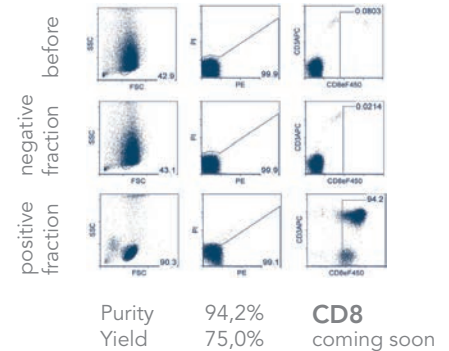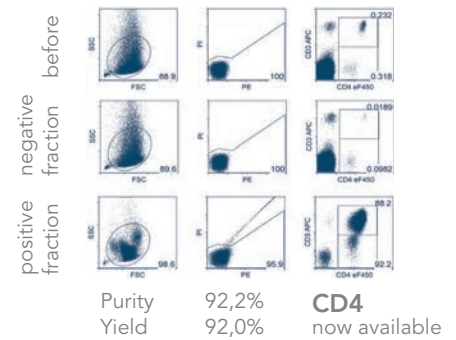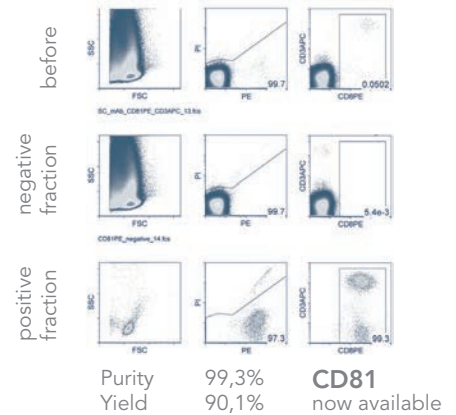

SELECTED CELL TYPES (E.G. PBMCs, T CELLS OR B CELLS) FROM WHOLE BLOOD

TECHNOLOGY


FABian® is equipped with a pre-made column filled with a Strep-Tactin® coated cellgrade agarose matrix and max. 11 tubes containing the required selection reagents.

FABian® starts the automated cell selection process by adding a specific Strep-tagged Fab fragment to the matrix. Subsequently, FABian® passes whole blood or other single-cell suspensions through the column. Target cells bind to the matrix based on the binding specificity of the loaded Fab fragment and non-target cells are washed away.


In the final step, FABian® adds biotin to the column, which causes elution of the Strep-tagged Fab from the matrix as well as dissociation of the Fab fragments from the cell surface.

The label-free authentic target cells are now selected in high yields and purity.


Remarkably, FABian® enables the selection of even small cell populations with high purity directly from whole blood.

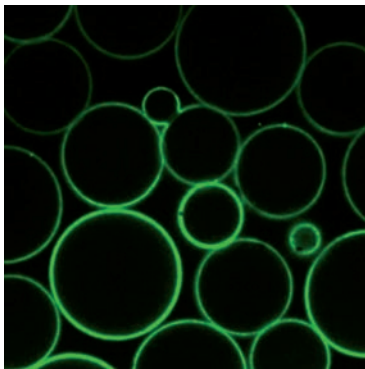

matrix loaded with fluorescently labeled Strep-tagged Fab fragments

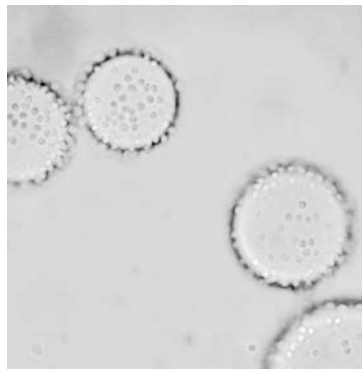

matrix loaded with target cells

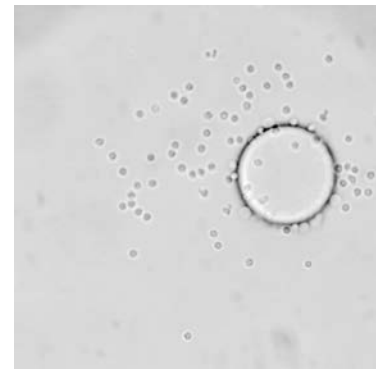

after adding biotin: target cells are released from the matrix

SERVICES

IBA GmbH continuously expands its cell selection applications and provides fully customized services for Fab fragment developments. If you require specific Fabs for your cell selection application that are not available yet please contact us.

www.fabian-online.com

IBA GmbH Rudolf-Wissell-Str. 28 · 37079 Goettingen · Germany · Phone +49 551 50672-0 · Fax +49 551 50672-181 · info@iba-lifesciences.com

IBA Sales Office Leipzig Bosestraße 4 · 04109 Leipzig · Germany · Phone +49 341 993898-29 · Fax +49 341 993898-19 · info@iba-lifesciences.com